

CRISPR Screening for Target Discovery in Primary Human Monocyte-derived Cells Using the VISTA Platform


Daniela Flehberger, Sofia Buianova, Emma Langguth, Johanna Irnstorfer, Stefanie Fellingner, Henrik Schmidt, Adam Krejci, Tilmann Buerckstuemmer, Anke Loregger, Lukas Badertscher, and Sumit Pawar

Myllia Biotechnology GmbH | Vienna, Austria


About Myllia's VISTA™ platform

Macrophages and dendritic cells are central players in the innate immune system. Their dysregulation is implicated in a wide range of diseases, including autoimmunity, inflammation, cancer, and infectious diseases. However, novel drug target identification in these cells has been difficult because these cells were - until now - refractory to high-content CRISPR screening approaches. At Myllia, we have solved this problem and have developed the **Versatile Integrated Screening for Target Analysis - VISTA™** platform, enabling CRISPR screens in primary human myeloid cells. The platform links CRISPR perturbation to transcriptomic responses, thus providing high-content functional readouts of novel gene targets involved in regulating many processes, including the tumor-immune interface relevant for immuno-oncology. **Here, we present a pooled CRISPR screen to dissect IFN-γ signaling in monocyte-derived macrophages (MDMs) and dendritic cells (MDCs).** Our platform thus enables large-scale CRISPR screens with complex readouts to support drug discovery campaigns.


1 The VISTA™ platform

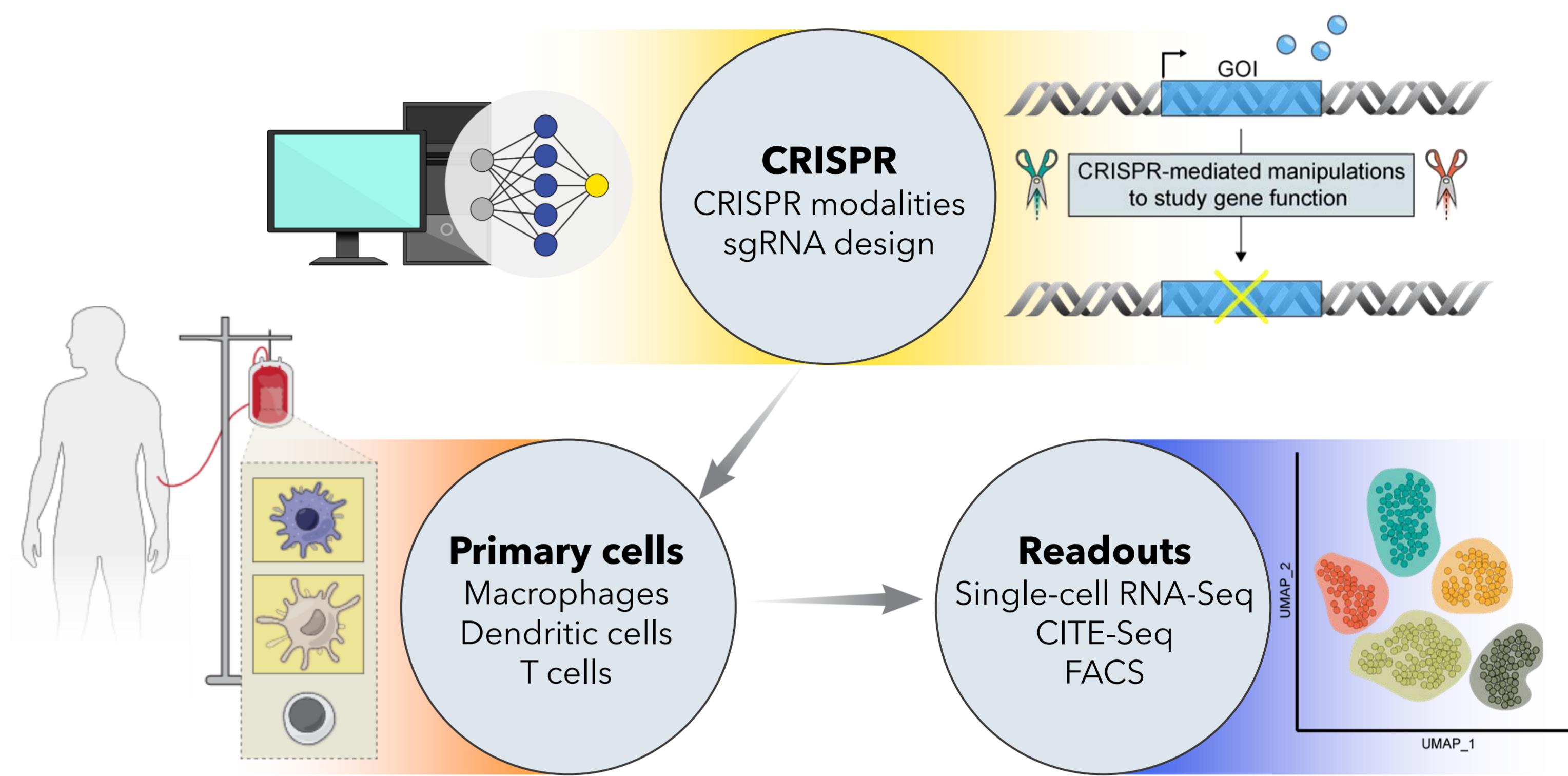


2 Macrophage polarization mechanisms

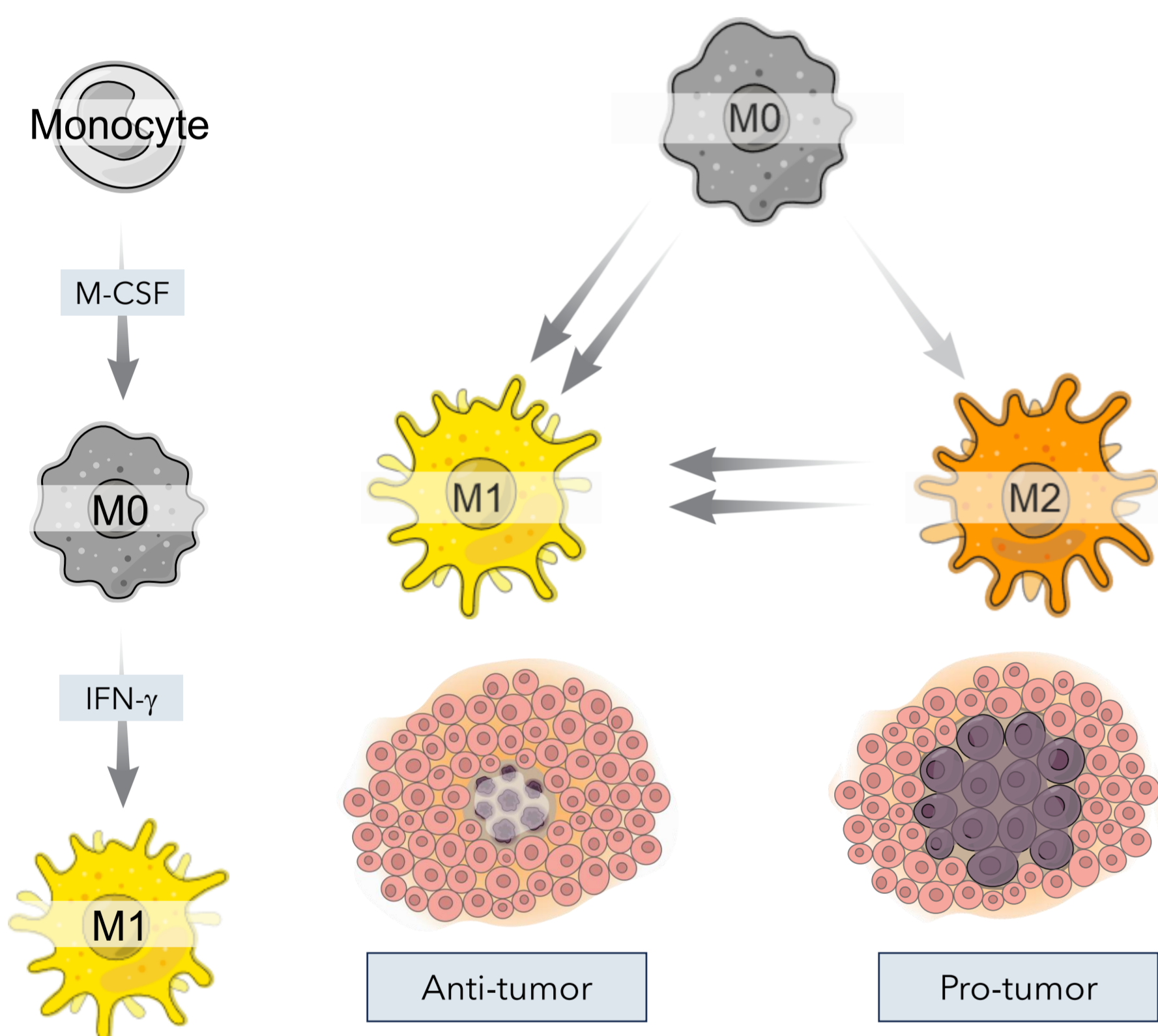


3 CRISPRi screens in macrophages

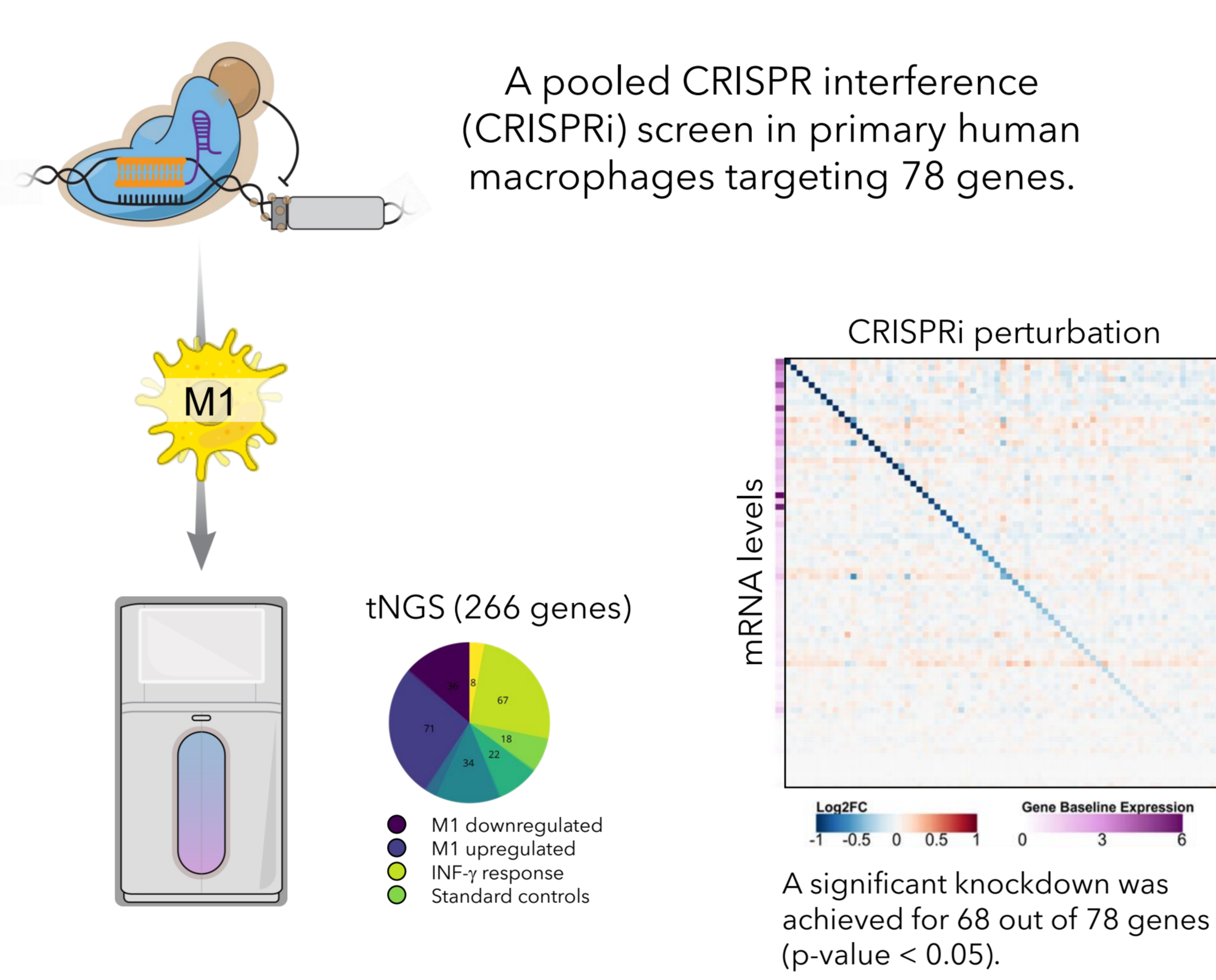


4 Polarization states resolved by scRNA-Seq

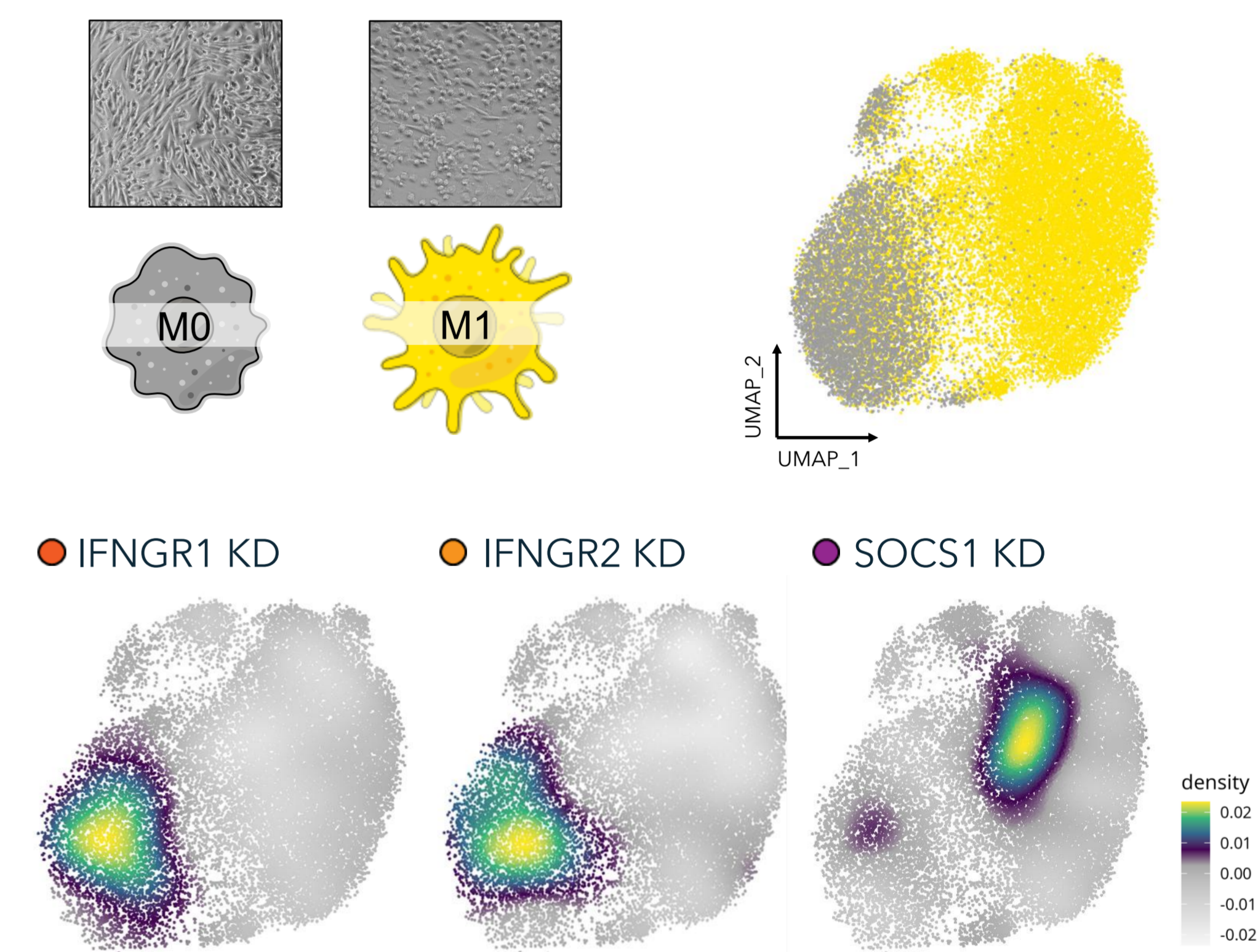

5 VISTA™ identifies key macrophage regulators and signatures

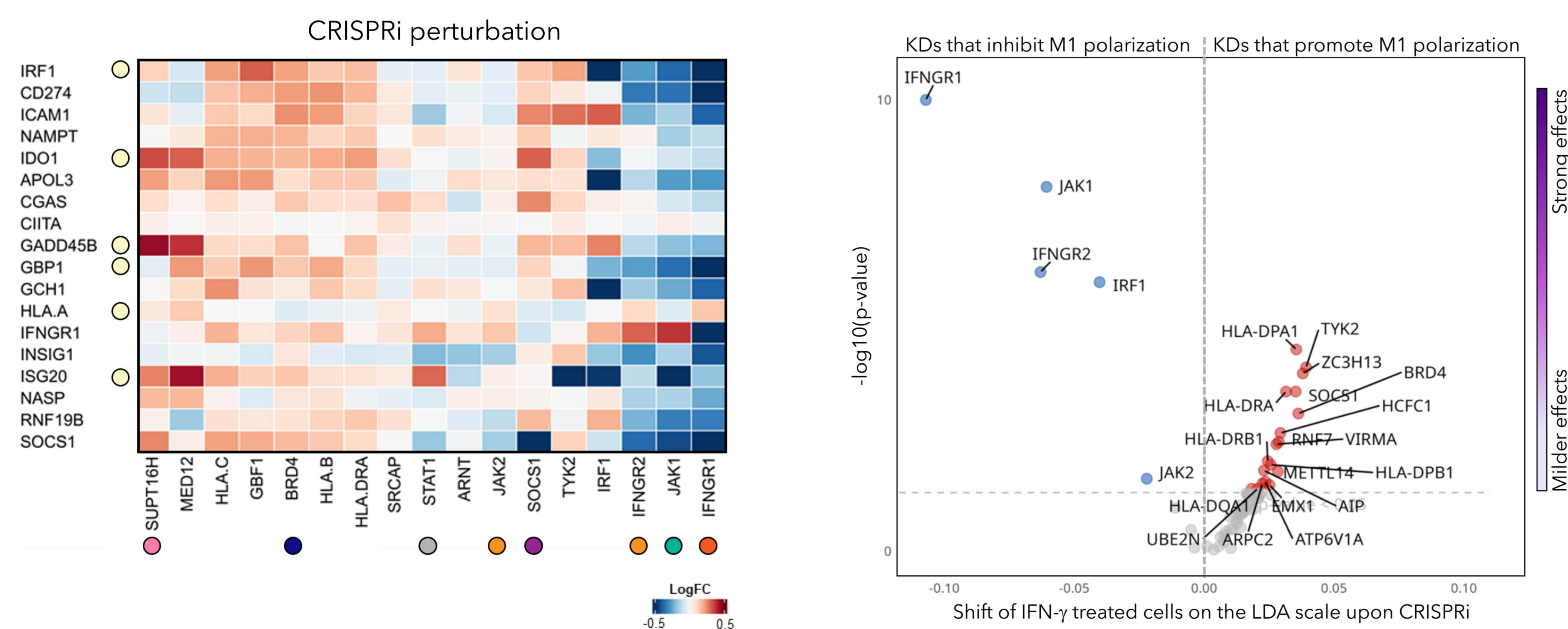

6 Perturbations affecting IFN-γ signaling

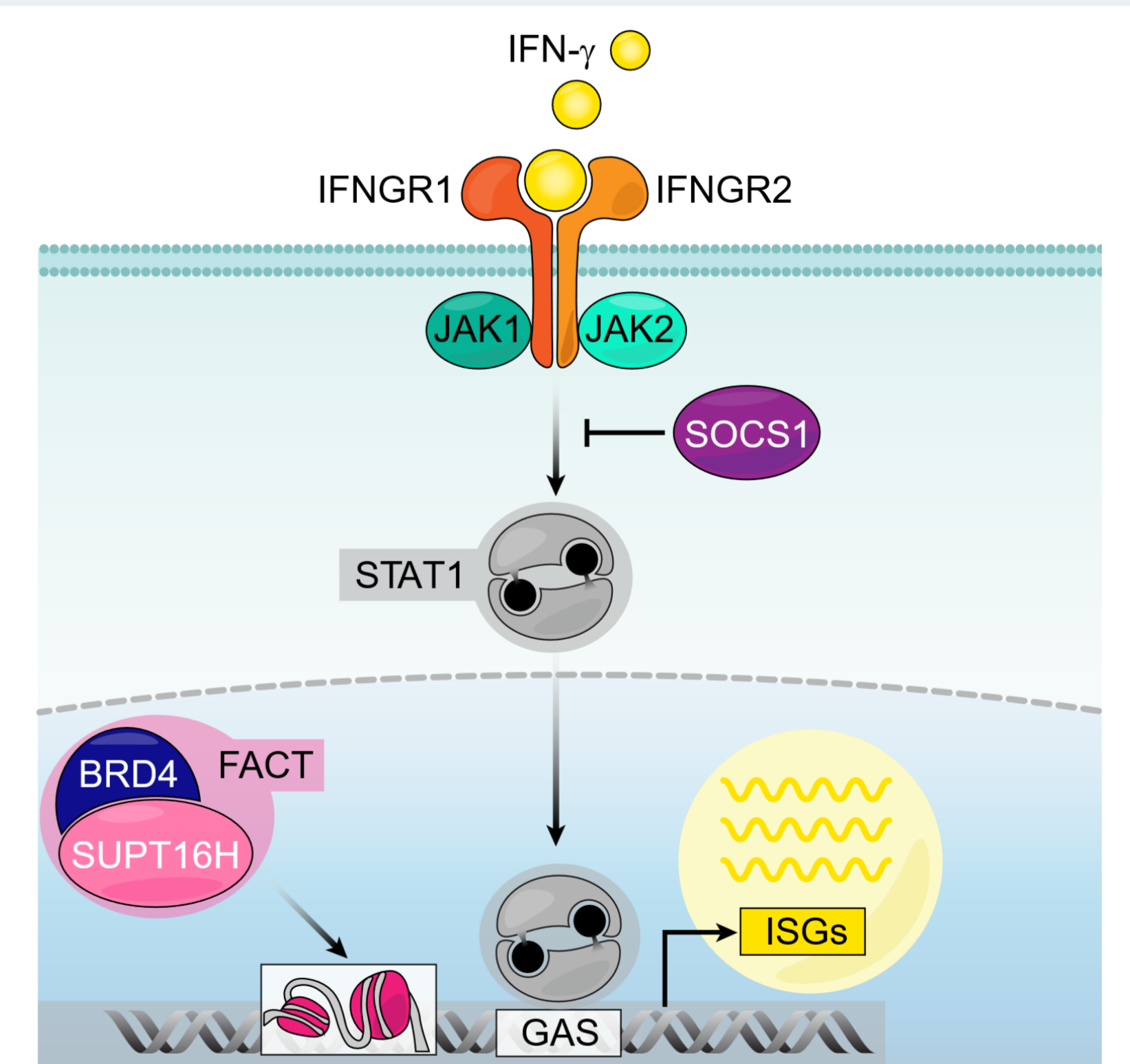


Conclusions and Outlook

- 5 positive and 20 negative regulators were identified that modulate M1 polarization via IFN-γ signaling
- The VISTA™ platform enables pooled CRISPR screens combined with single-cell RNA-Seq in various primary human cell types, including macrophages, dendritic cells (and T cells)
- Key features of Myllia's VISTA™ platform include:
 - Unbiased drug target discovery
 - Target identification in blood-derived primary human cells
 - Targets with higher chance of translating *in vivo*

We are looking for a strategic partner to initiate drug target discovery programs across several disease areas in primary human macrophages and dendritic cells, incl. autoimmune disease, inflammation and oncology.

Contact us

Myllia Biotechnology
www.myllia.com

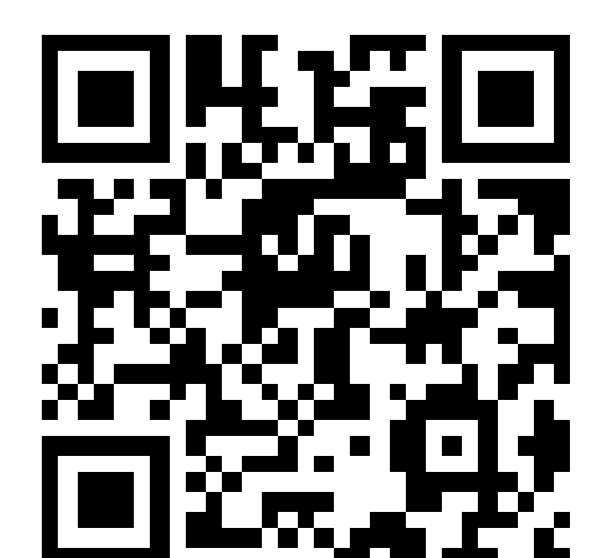

E: info@myllia.com

